


Astral™

Inoculation System

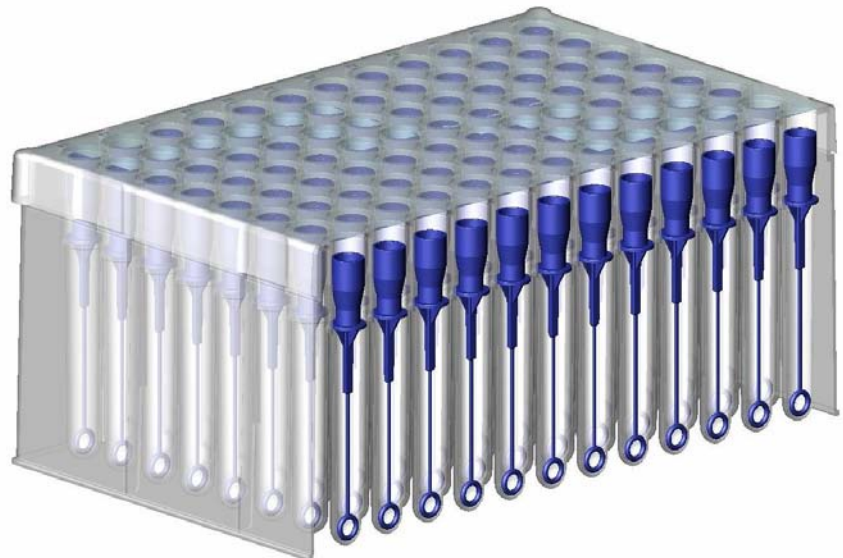

BIO PLAS, Inc.,
4340 Redwood Hwy, Suite A15, San Rafael, CA 94903 USA
Tel: (415) 668-8917 Fax: (415) 668-8354
e-mail: info@bioplas.com web: www.bioplas.com

Astral™ Inoculation System

The disposable BIO PLAS Astral Inoculation System is color coded and consists of the following; 1 uL yellow Astral Loop, 10 uL blue Astral Loop and white Astral Streaking Needle.


During manufacturing, these Astral Loops are inserted into racks which have 96 individual tubular chambers. When the 96 chambers are filled with the Astral Loops they are then hermetically sealed with a thin plastic membrane. A dust cover is placed over this 96 chamber rack, heat shrunk and sterilized by irradiation.

Additionally all Astral Loops are certified to be free of RNase, DNase, DNA, pyrogens and PCR inhibitors.

Every case of 10 racks of 96 Loops is supplied with a free Astral Handle.

The handle is designed to be used with one hand and will easily puncture the membrane attaching itself to the hub of the Astral Loop. It may then be withdrawn from its individual sterile chamber ready to be used for inoculation or streaking. When the inoculation or streaking process is complete, the Astral Loop or Astral Streaking Needle may then be ejected off the Astral Handle.

This easy to use and ruggedly designed Astral Handle may be autoclaved repeatedly, without damage to the handle.

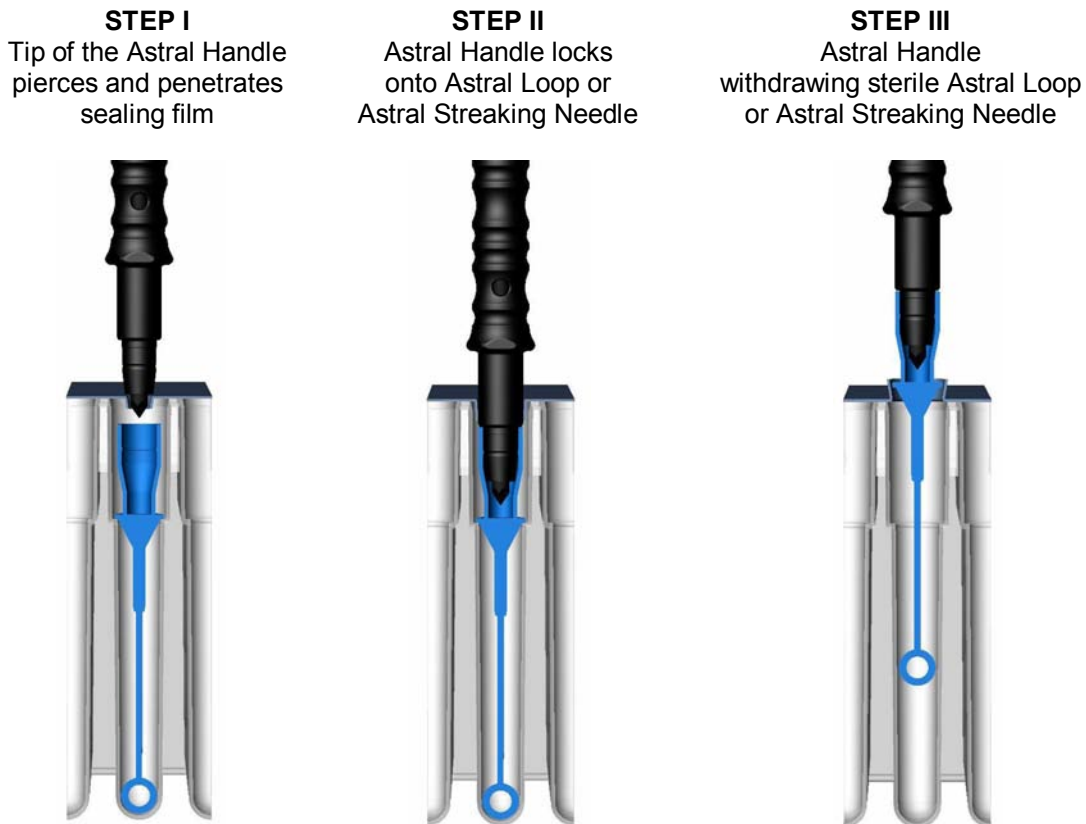


This illustration shows the three easy and rapid steps used to extract a sterilized 10uL Astral Loop from its sealed chamber.

BIO PLAS Inc.,

Phone: (415) 472-3777 Fax: (415) 472-3758

email: info@bioplas.com website: www.bioplas.com

Procedures: All Astral Loops may be used in quantitative procedures such as sampling, urine counts, serial dilutions, and bacteriological inoculation.

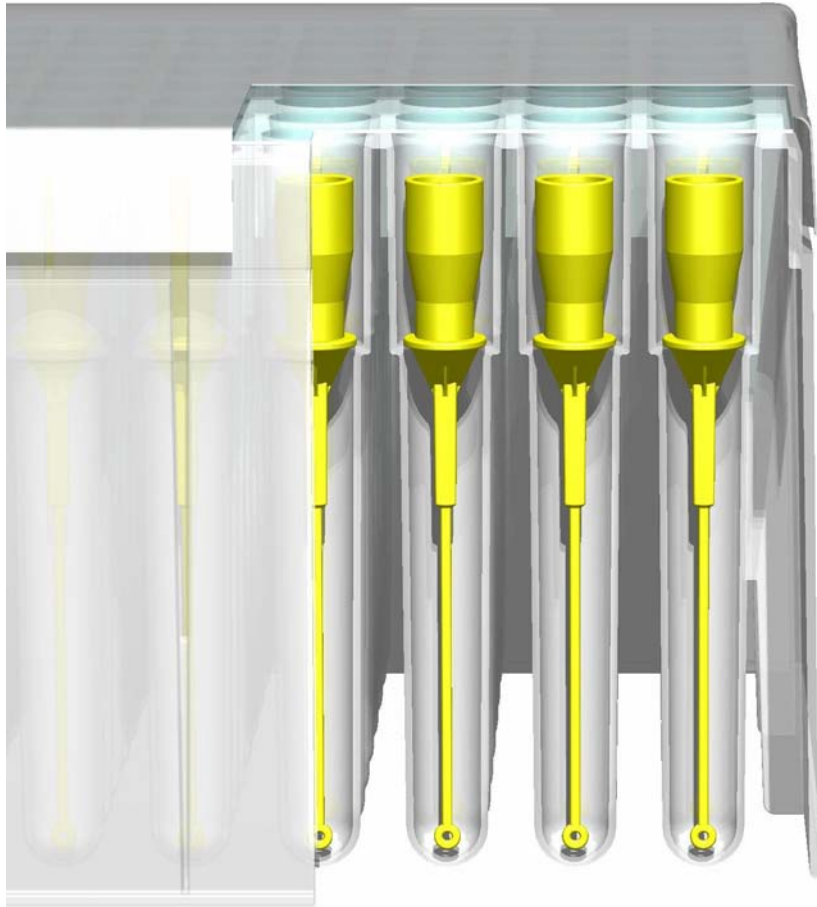

**1uL Astral
Inoculating Loop**
Catalog Number 7000

SAFETY: Astral Loops eliminate the metal or wire loop flaming step and the hazards of spattering, cross-contamination, and the risk of infectious pathogens caused as a result of aerosols, and the hazards of open flame.


CONVENIENCE: All Astral Loops may be quickly extracted from the sterilized rack by using the Astral Handle. These Astral Loops are flexible enough to facilitate streaking but rigid enough to penetrate the media for inoculation.

TIME SAVINGS: The Astral Inoculation System allows the technologist to inoculate prepared plates at an extremely rapid rate in any laboratory setting or condition, i.e., anaerobic chambers, fume hoods, satellite labs or, in the field.

ECONOMY: The savings can be substantial; fewer plates used, guaranteed sterility and cleanliness, zero cross-contamination and speed of use.

BIO PLAS, inc.

Astral™ Inoculation System

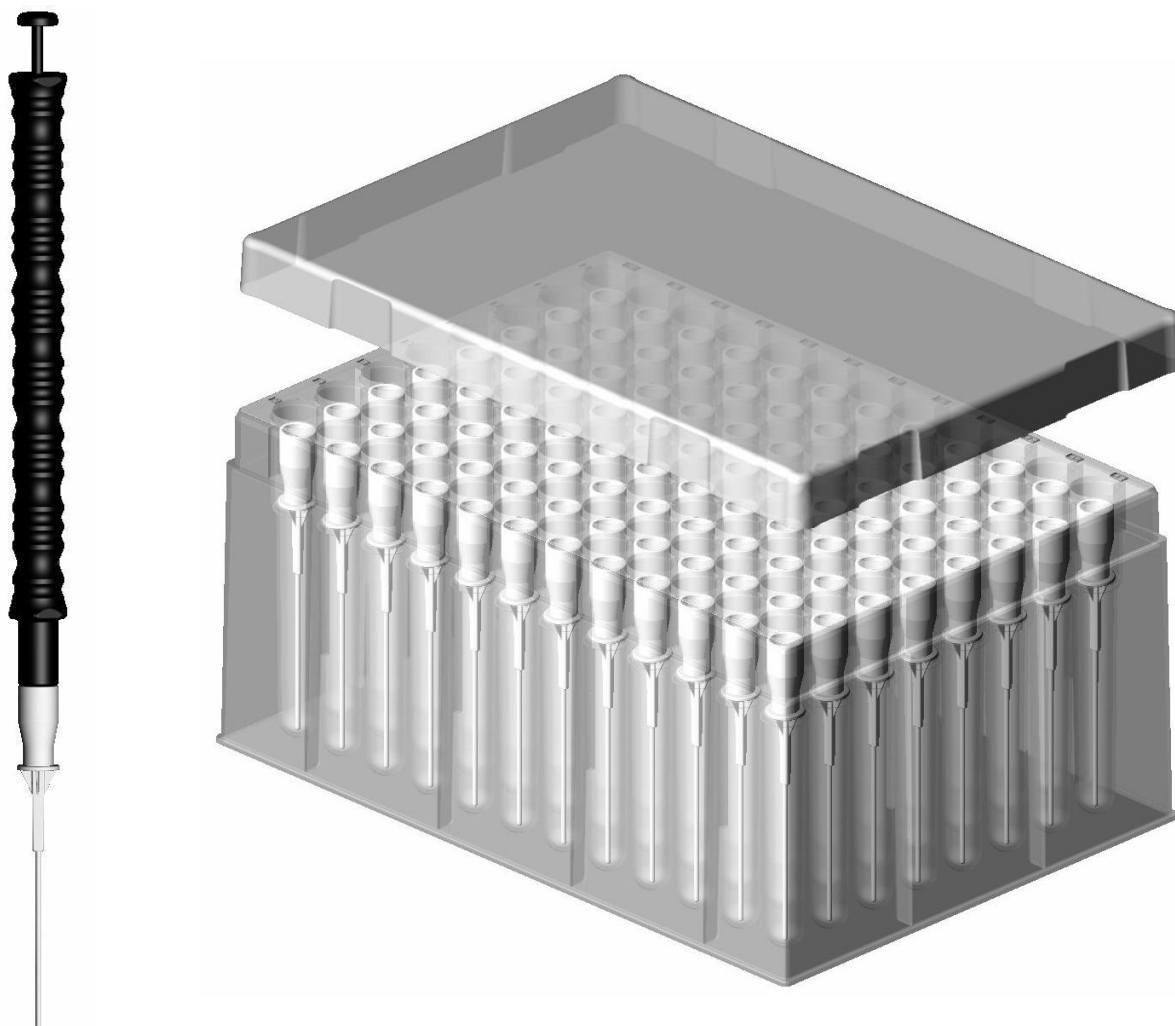

Ordering Information

Catalog #	Description	Pkg. Qty.	Pkgs./ Case
7000	1uL Astral Inoculation Loop, Yellow	96/Rack	10
7010	10uL Astral Inoculating Loop, Blue	96/Rack	10
7020	Astral Streaking Needle, White	96/Rack	10
7050	Astral Handle	1	1
<i>Astral™ is a Trademark of BIO Plas, inc.</i>		Minimum Order One Case.	
<i>*Patent Number: 5,330,899, Patent Pending</i>			

BIO PLAS Inc.,

Phone: (415) 472-3777 Fax: (415) 472-3758

email: info@bioplas.com website: www.bioplas.com